V CARE SteriControl

DEUTSCH

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ENGLISH

TRYPTIC SOY BROTH (TSB) MEDIUM FOR A CONTROLLED WORKING ENVIRONMENT

INFORMATION

- Growth promotion tested
- Sterility confirmed
- Shelf life 12 months from date of production
- 20 ml glass and PET(G) bottles

- V CARE SteriControl enables an easy control of the aseptic working environment and medium handling. The product enables an unmatched level of control over the sterile liquid handling processes used in ART and can be used alone or in addition to sedimentation plates.
- V CARE SteriControl is handled like a culture medium in the working area, thereafter incubated and visually controlled for turbidity after two weeks.
- V CARE SteriControl ensures the traceability of sterile working conditions.

WARNING

- Do not use after expiry date.
- Do not re-sterilize after opening.
- VITROMED recommends only handling of media with aseptic techniques (e. g. LAF-bench ISO Class 5).
- Do not use the product if any signs of microbial contamination are noticed.
- Do not use if discolored, turbid or a precipitate is observed.
- Do not use if seal of the container is damaged upon delivery.
- Do not use medium that shows any evidence of particulate matter, cloudiness, or color change.
- Never place biological materials, such as sperm, oocytes or embryos in V CARE SteriControl!

PRECAUTIONS

- Only for intended use.
- Leave product in its original packaging.
- Discard minimal amounts of excess medium remaining in the bottle.

INSTRUCTIONS

WORKPLACE CONTAMINATION PROTOCOL

Use V CARE SteriControl as indicator of workplace sterility or bioburden (the number of microbes present):

1. Open the bottle in the working area.
2. Keep the bottle open while the work is being performed.

Caution: Never place biological materials, such as sperm, oocytes or embryos in V CARE SteriControl!

3. Close the bottle.
4. Incubate the bottle at 20-35 °C for 2 weeks. (Ideally, the first week at about 22 °C and the second week at about 33 °C).
5. Evaluate the media in the bottle visually: for sterility after 2 weeks; for bioburden daily.

HANDLING AND PLASTICWARE CONTROL PROTOCOL

Handle V CARE SteriControl like a culture medium, but without biological materials in the working area:

1. Open the bottle in the working area.
2. Conduct all your regular work steps without biological material and use V CARE SteriControl instead of your regular ART medium.

Caution: Never place biological materials, such as sperm, oocytes or embryos in V CARE SteriControl!

3. Keep the bottle open while the work is being performed.
4. Close the bottle.
5. Incubate the bottle at 20-35 °C for 2 weeks. (Ideally, the first week at about 22 °C and the second week at about 33 °C).
6. Evaluate the media in the bottle visually: for sterility after 2 weeks; for bioburden daily.

CONTROL OF MEDIA CONTAMINATION PROTOCOL

Use V CARE SteriControl to evaluate contamination of culture and handling media

1. Open the bottle in the sterile working area.
2. Quickly add up to 1,0 ml of the medium under investigation into the bottle of V CARE SteriControl!

Caution: Never place biological materials, such as sperm, oocytes or embryos in V CARE SteriControl!

3. Close the bottle.
4. Incubate the bottle at 20-35 °C for 2 weeks. (Ideally, the first week at about 22 °C and the second week at about 33 °C).
5. Evaluate the media in the bottle visually: for sterility after 2 weeks; for bioburden daily.

PRODUCT SPECIFICATIONS

- Sterility: sterile (no growth)
- Growth promotion test: passed
- Visual control: passed

STERILE A Sterilized using aseptic processing techniques

EVALUATION OF THE RESULTS

The media appears turbid: contamination.

The media appears clear: no contamination.

Early growth indicates a high bioburden. The cleanliness of the workplace should be improved.

No growth after 14 days indicates a sterile process.

Growth can also be evaluated with the Tyndall effect using a Laser Pointer. The Tyndall effect is light scattering by particles in a colloid solution or in a very fine suspension. When the sample is not contaminated, no laser beam as light scattering can be seen. When there is contamination a light scattering by the fine particles is obvious.

V CARE SteriControl can be also used to do plating and sub-cultivation.

STORAGE INSTRUCTIONS AND STABILITY

- Do not use after expiry date.
- Store between 2 °C to 37 °C.
- The product can be used for a period of up to 7 days after opening, provided that it is handled under aseptic conditions.
- Keep away from sunlight.

FEEDBACK AND OBLIGATION TO INFORM

Thank you for purchasing our product. If you have any suggestions for us to improve our products, please send them to qc@vitromed.com.

Report any serious incidents that have occurred with the product immediately by email to qc@vitromed.com and the competent authority of the Member State in which the user and/or patient is established or treated.

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Technischer Support **VITROMED GmbH**